

# PERVATECH<sup>®</sup>

## Datasheet

# FLAT SHEET POMS MEMBRANE

Flat sheet POMS (polyoctylmethyl siloxane) membranes have organophilic (hydrophobic) characteristics, which means that the organic constituent of the feed passes preferentially through the membrane.

### MEMBRANE ELEMENT

Geometry	Flat sheet
Dimensions	297 x 210 mm per flat sheet (package of 3 pcs.)
Substrate material	PET
Intermediate layer	Type of PI
Top layer	POMS

### OPERATIONAL WINDOW

Temperature	Maximum 70 °C (short-term 80 °C)
Pressure	Maximum 5 bar
pH	3-10
Pre-filtration	10 µm cartridge filter

### HANDLING, CLEANING AND STORAGE

#### **Handling**

- Wear clean gloves to prevent contamination with fungi.

#### **Cleaning**

- Clean the element by flushing with water to which a non-ionic detergent (10 ppm KOH) has been added. Alternatively, enzymatic solutions may be used, depending on the feed composition.
- In case of food & additives processing, please contact us for alternatives.

## Sterilization options

Steam

106-108 °C

Flushing

With ethylene oxide or 100 % ethanol

## Storage

Storage before use

Out of direct sunlight, at room temperature, < 70 % relative humidity

Storage after use (short-term)

In a mixture of water and 10-15 % IPA or in a solution of water with 2500 ppm sodium metabisulfite

Storage after use (long-term)


In a solution of water with 0.7 % benzalkonium

## EXAMPLES OF APPLICATIONS WITH FLAT SHEET ORGANOPHILIC (HYDROPHOBIC) POMS MEMBRANES

- Recovery or extraction of organics (in biotech or biotech-related food applications) from natural feed streams like fruit juices, wine, beer, coconut oil, essential oils (carvone, limonene) and in combination with fermentation
- Removal of ethyl alcohol (and other alcohols) from wine and beer
- Upgrading of reverse osmosis permeate in juice production
- Combination with bioreactor in production of alcohols (ethanol, IPA, butanol), ABE (acetone, butanol, ethanol), aldehydes, flavor production as well as acid production
- Removal of volatile organic compounds (VOC)

In general, POMS membranes are more open than PDMS membranes. However, in some solutions, especially  $\gg$  C2 molecules and low concentrations, POMS performs better than PDMS. This especially holds for phenolic types of molecules.

Typical example of pervaporation (PV) process in combination with bioreactor:

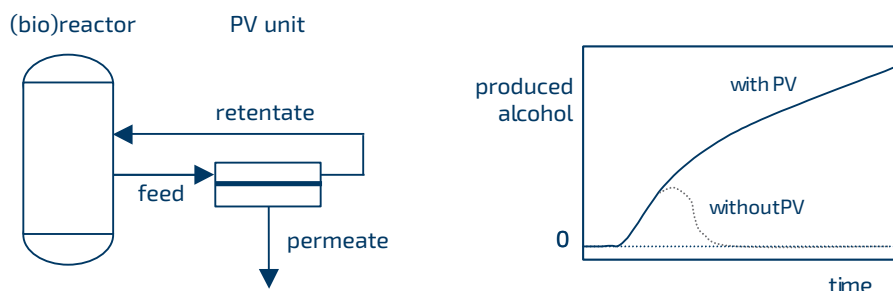

The image shows the principle of a membrane reactor for continuous product recovery (alcohols, aromas). The system with pervaporation continuously produces alcohol while other systems stop when inhibiting amounts of alcohol (or another inhibitor) have been reached. In some cases, a pretreatment is necessary.

Heliumstraat 11  
7463 PL Rijssen  
The Netherlands  
T +31 548 530 360  
info@pervatech.nl  
pervatech.com